



PickPen[®] 8-M

the unique magnetic tool for parallel sample processing

Rapid isolation of nucleic acids

Rapid purification and fractionation of proteins

Use with optimized QuickPick™ kits and ready-to-use DNase/RNase free tips



FAST

- move particles, not liquids to reduce number of steps
- process samples using 96-well plate format

EFFICIENT

- simultaneous processing of 8 samples
- eliminate need for shared equipment
- increase throughput without major investment

EASY

- simple to learn, operate and maintain
- hand-held, carry PickPen[®] with you
- excellent usability through unique ergonomic design

GENTLE

- ensure high sample quality through rapid handling
- avoid centrifugations and vacuum manifolds



PickPen® 8-M: principle of operation

The key to PickPen® 8-M is its innovative magnet construction* and the elastic tips which fit closely over the magnets to minimize dead volume. This allows work with even the smallest sample volumes.

When locked in the extended position the magnets collect magnetic particles from 8 samples simultaneously. When the magnets are withdrawn, the particles are released into the next solution. This action can be repeated for as many steps as needed, using the same tips. Sample cross-contamination is eliminated as tips are always changed between samples.



The proprietary PickPen® tips are delivered in a unique ready-to-use pack with 8 DNA/RNase-free tips. All QuickPick kits contain tips.

PickPen® 8-M: efficient and cost effective

PickPen® 8-M is designed for higher throughput and 96-well plate format. Increase efficiency without the investment needed for dedicated systems while maintaining all the benefits of the PickPen® technology. The proprietary ergonomic design makes the tool comfortable and easy to use. PickPen® is optimized for use with Bio-Nobile QuickPick kits, but can be used for other magnetic particle based applications as well.



PickPen® products:		BN QuickPick™ kits:	
PickPen® 8-M	23008	mRNA purification kits	
PickPen® tips 8-pack, DNase/RNase free	34008	DNA purification kits	
PickPen® 96 tips	34096	Protein purification kits	
PickPen® Tip box (no tips)	34051	Protein fractionation kits	
PickPen® 1-M	23001		

Distributed by:



info@bio-nobile.com
Tel. +358-2-410 1100
P.O. Box 36, FIN-20521 Turku, Finland

PickPen is a registered trademark and QuickPick a trademark of Bio-Nobile Oy.

*patent pending